LabVIEW Anwender Treffen

Was ist neu in LabVIEW 2014

Marcel Krenz

<http://www.fhi-berlin.mpg.de/pc/labview>

06.11.2014

Fritz-Haber Institute, Berlin


Accurate Data Acquisition With
the Latest Technology

Extended Deployment Options
for Built-In Analysis Routines

13 Coding Productivity
Enhancing Features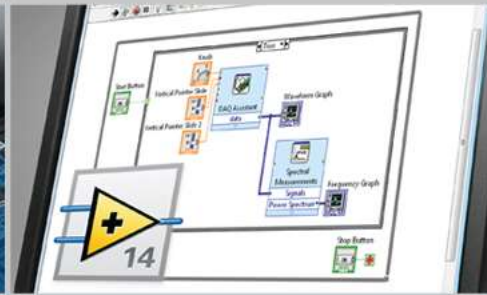


LabVIEW™ 2014

Acquire, analyze, and visualize any data set to make
informed decisions fast.

Expanded Interface Development
and Deployment Options

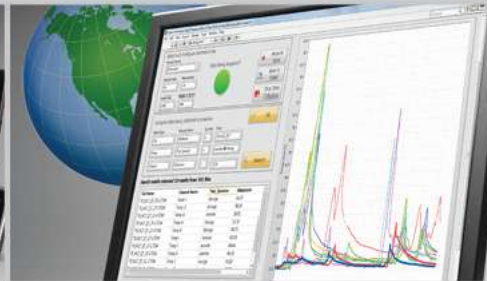
Global Solutions to Mine
and Search Data

Innovate With a
Community


Unrivaled Integration With the Latest Technology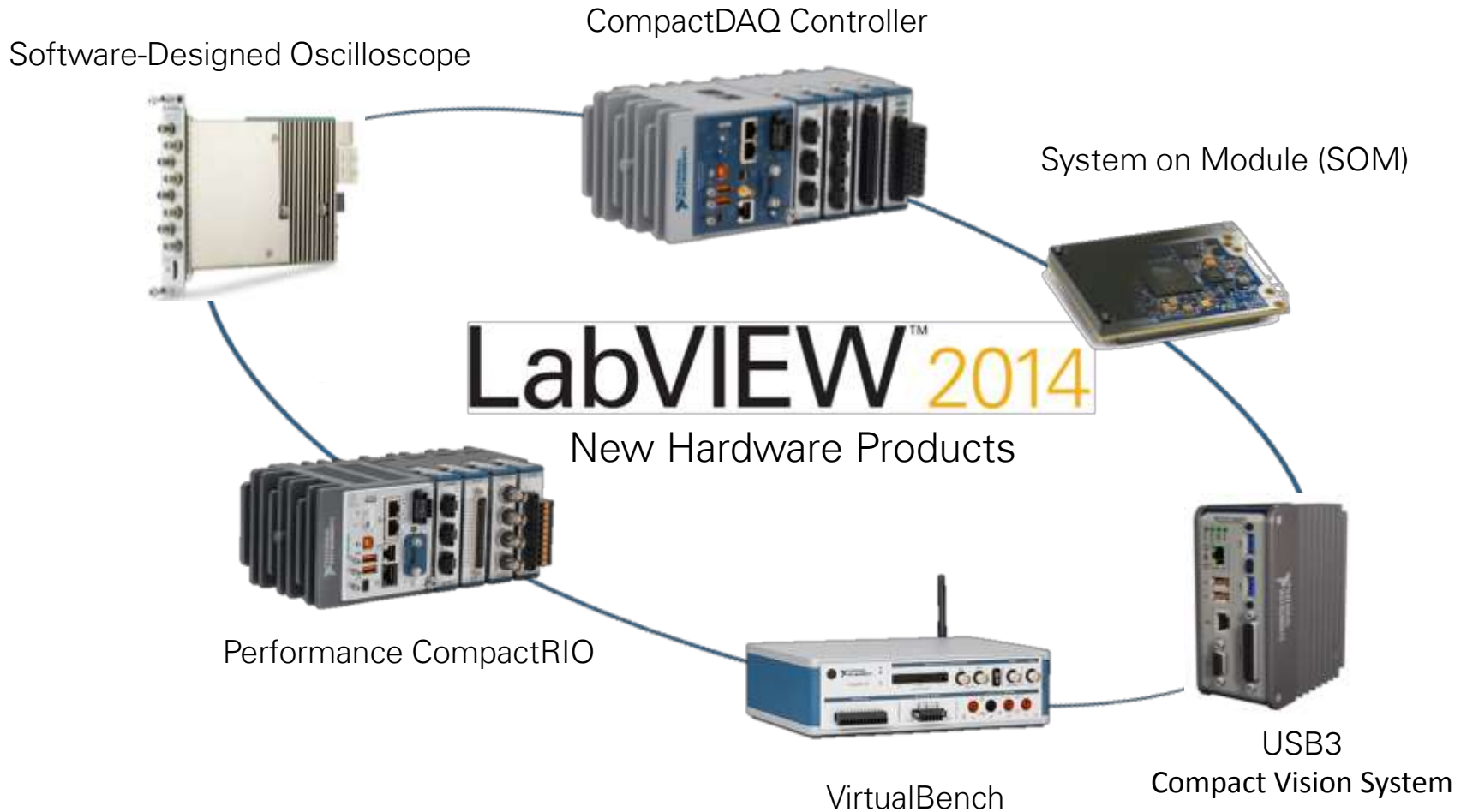

Simplified Connection to Real-Time Targets

Connect With USB

New NI Linux Real-Time targets use **Ethernet over USB** to connect to the host computer

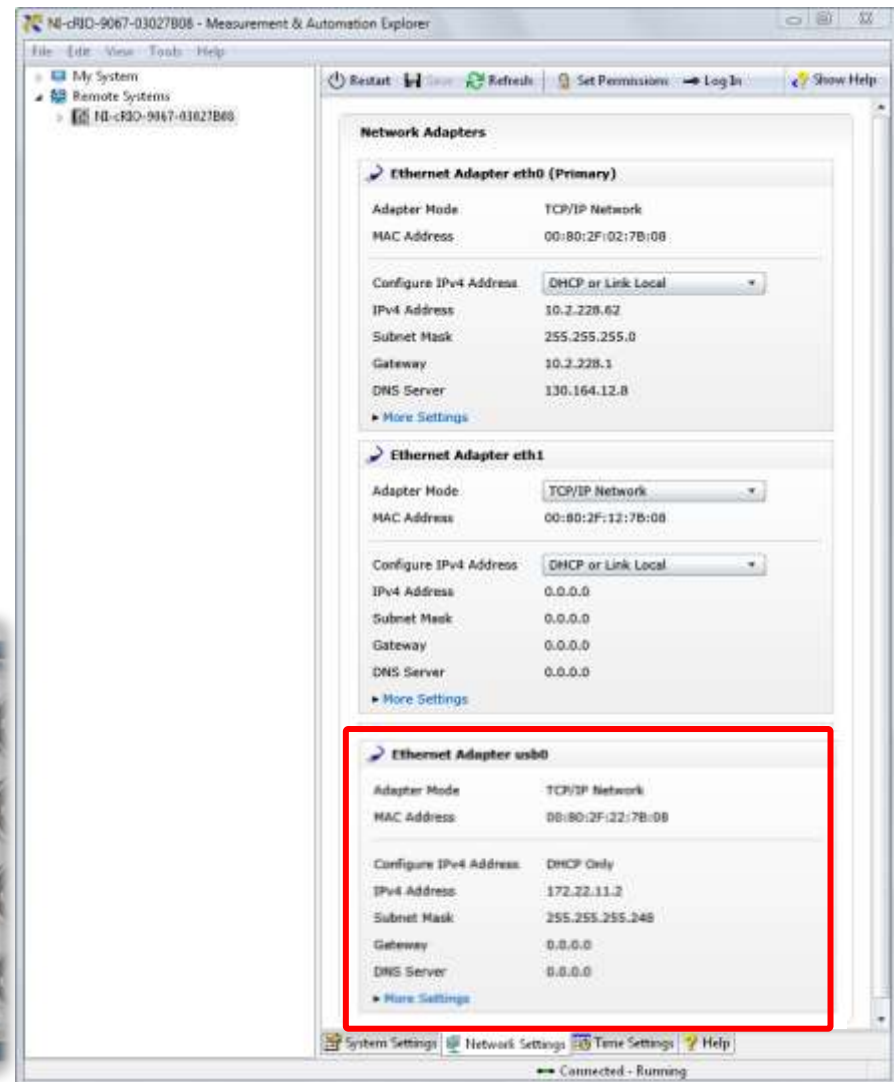Value Added to Existing LabVIEW Editions

Full Edition


Make advanced control applications through included **PID and Fuzzy Logic** palettes

Professional Edition

Trace low-level bugs, validate coding practices, and manipulate data through the now included:

- **VI Analyzer Toolkit**
- **Unit Test Framework Toolkit**
- **Desktop Execution Trace Toolkit**
- **Database Connectivity Toolkit**
- **Report Generation Toolkit**

ni.com

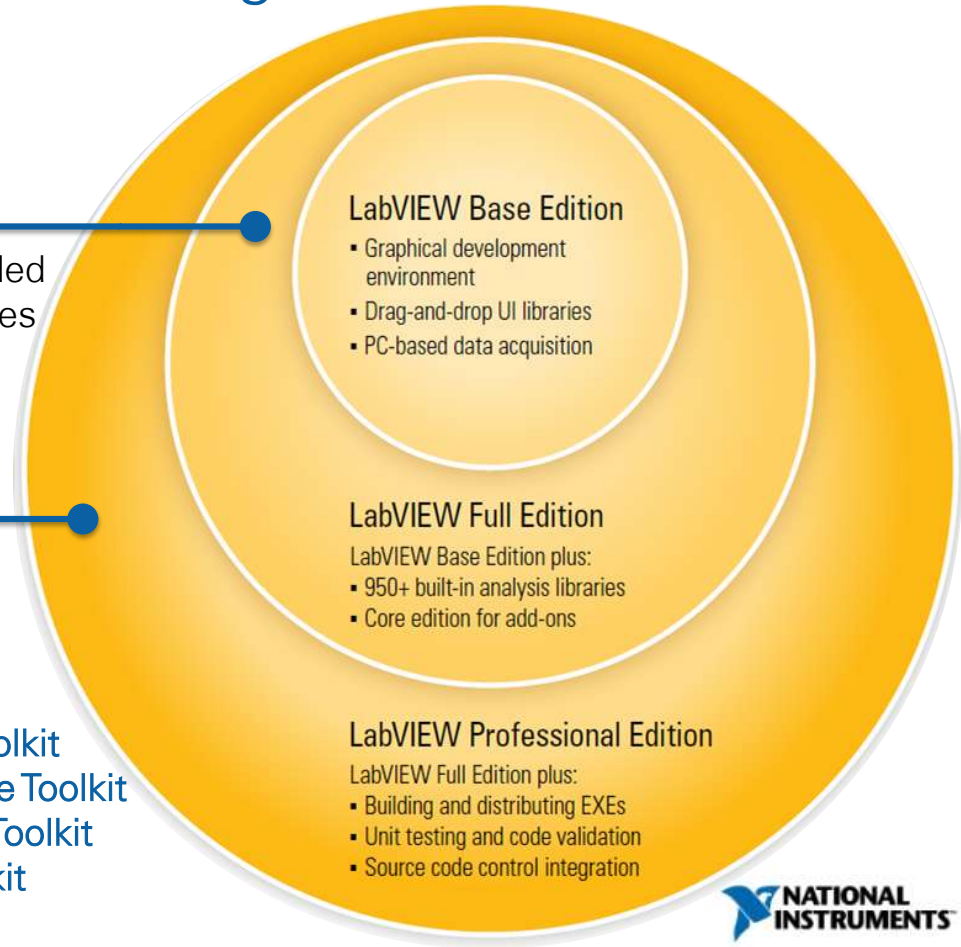New Taskbar Icon

- Easily differentiate between LabVIEW versions
- Suggested by 500+ [Idea Exchange](#) kudos

New Taskbar Icon

32-bit

64-bit

LabVIEW™ 2014

What's New in LabVIEW 2014

Kellis Garrett Wiseman

NI Field Sales Engineer

Troubleshoot Faster With Guided Error Debugging


Embedded resource links

Why do I get Error 43 when using the Open/Create/Replace File VI?


Primary Software: LabVIEW Development Systems>>LabVIEW Professional Development System
Primary Software Version: 8.2
Primary Software Fixed Version: N/A
Secondary Software: LabVIEW Development Systems>>LabVIEW Full Development System

Problem:
When I use the Open/Create/Replace File VI, I have the option to select a file path or cancel the operation. If I choose **Cancel** I get the error, **Error 43 occurred at Open/Create/Replace File in [VI Name]** .

1. Error Dialog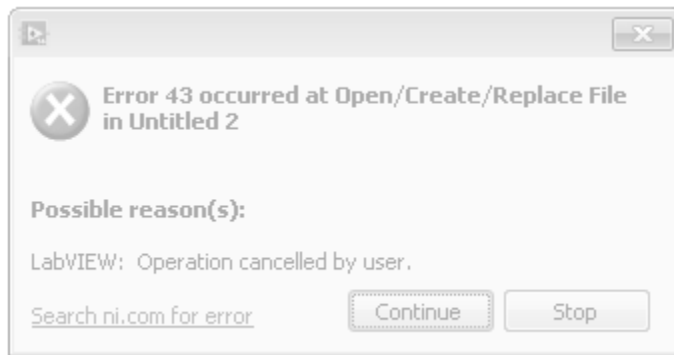

2. Search Result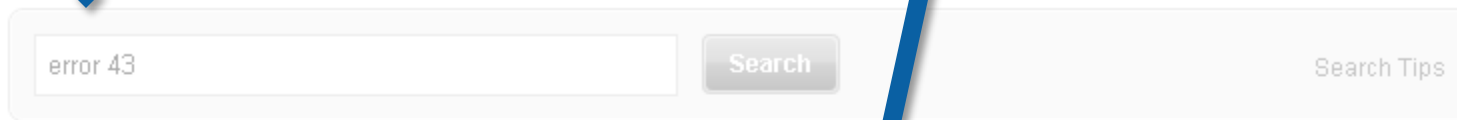

1-10 of 3,564 Show 10 20 30 results per page


1 2 3 4 5 6 7 8 9 10 Next

Why do I get Error 43 when using the Open/Create/Replace File.vi?

I get Error 43 when using the Open/Create/Replace File.vi? Primary Software: LabVIEW Development Systems>>LabVIEW... following error, ...

digital.ni.com/public.nsf/allkb/E78D4D60B08E4CCE862572970058B706

3. Solution

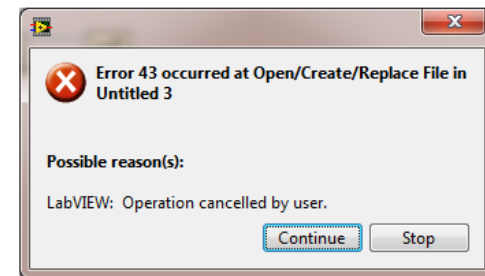
Time-Saving Development Tools

Automatic shift register creation

LabVIEW Idea Exchange

- For Loops output arrays by default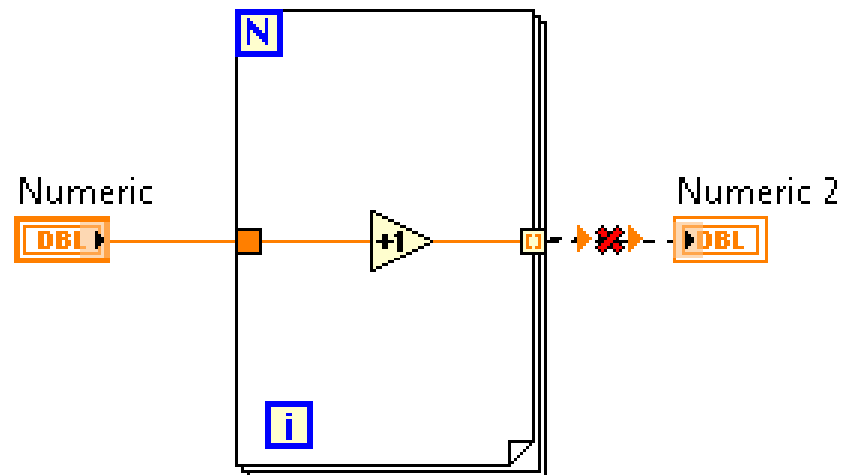

LabVIEW 2013


Time-Saving Development Tools

Automatic shift register creation

LabVIEW Idea Exchange

- For Loops output arrays by default
- Context-sensitive to wired outputs
- Shift register creation matches logical flow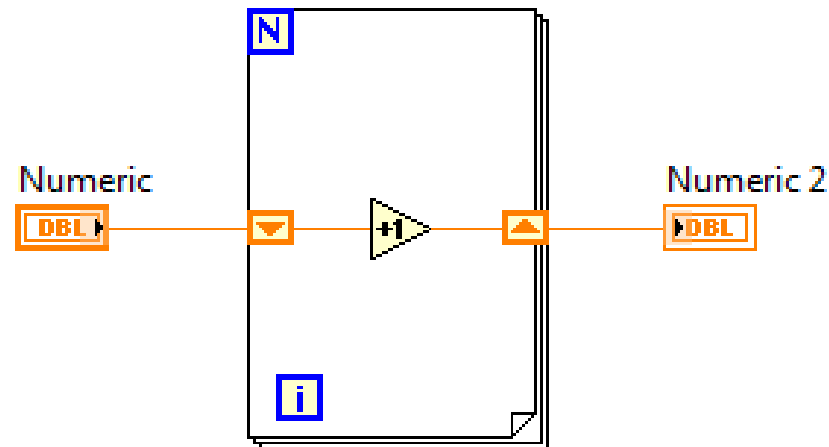

LabVIEW 2014


Time-Saving Development Tools

Replace case structure tunnel with selector

LabVIEW Idea Exchange

LabVIEW 2014

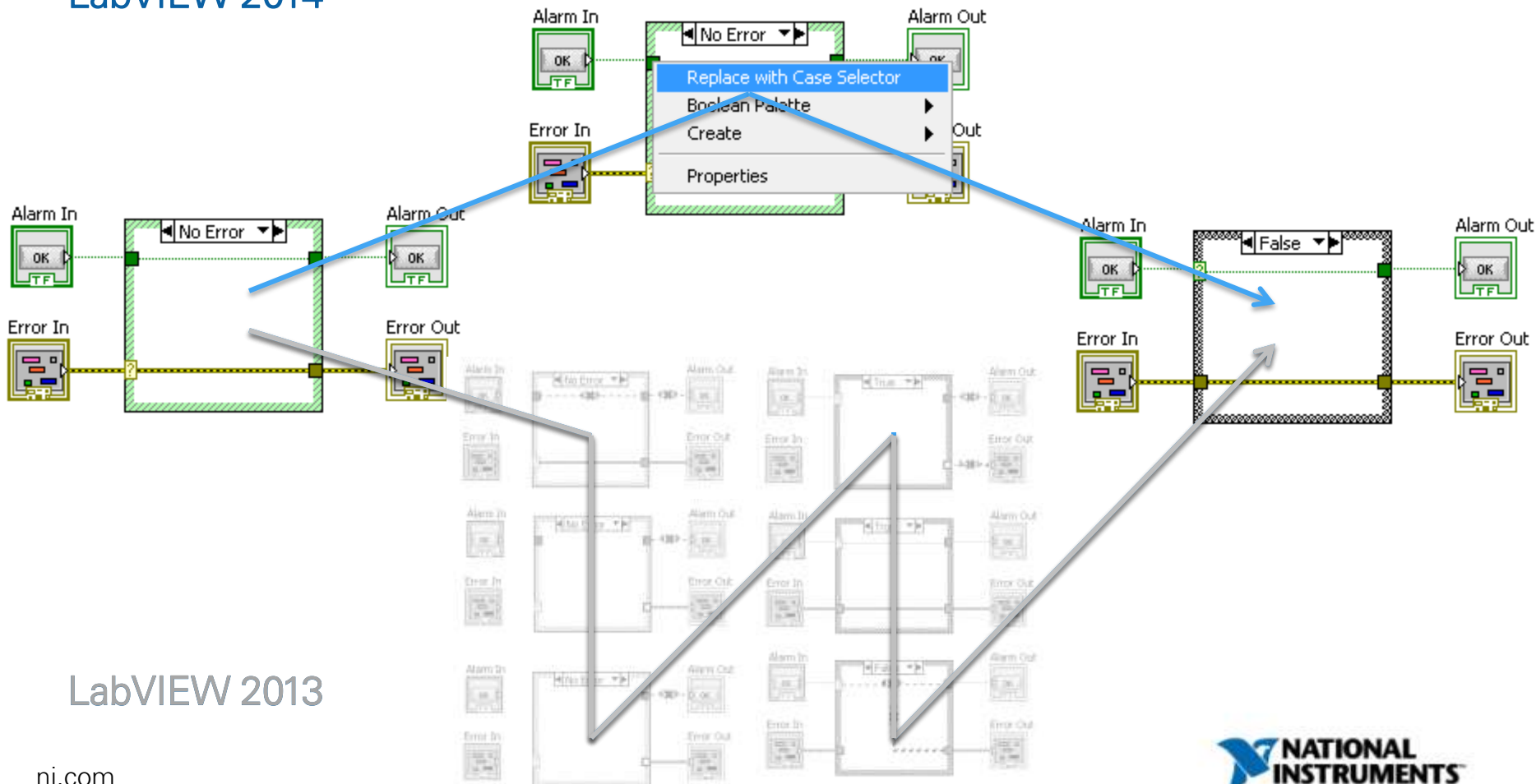
LabVIEW 2013

Time-Saving Development Tools

Auto-Wire Quick Drop shortcut

LabVIEW Idea Exchange

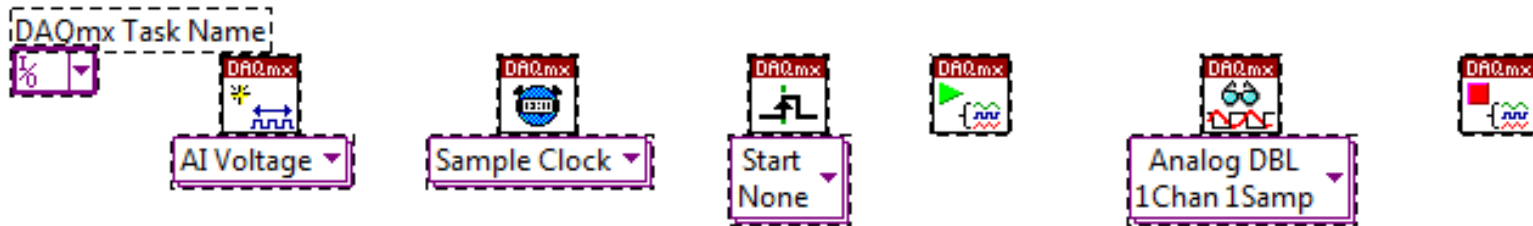<Ctrl-Space>
<Ctrl-W>

Data Dashboard for LabVIEW 2.2

Alternate Servers

Now you can set up one dashboard to monitor multiple targets and use the drop-down menu to switch where the data is streaming from while the dashboard is running

Multiplot

Data Dashboard now supports 2D array of numerics as a data type for charts and graphs

Android Tablets



Data Dashboard on Android tablets now supports monitoring and control applications and free canvas customizations

Transparency

Adjust the transparency of images, controls, and indicators on your dashboard

LabVIEW 2014 Real-Time With Embedded UI

Simplify system complexity by implementing a local HMI on CompactRIO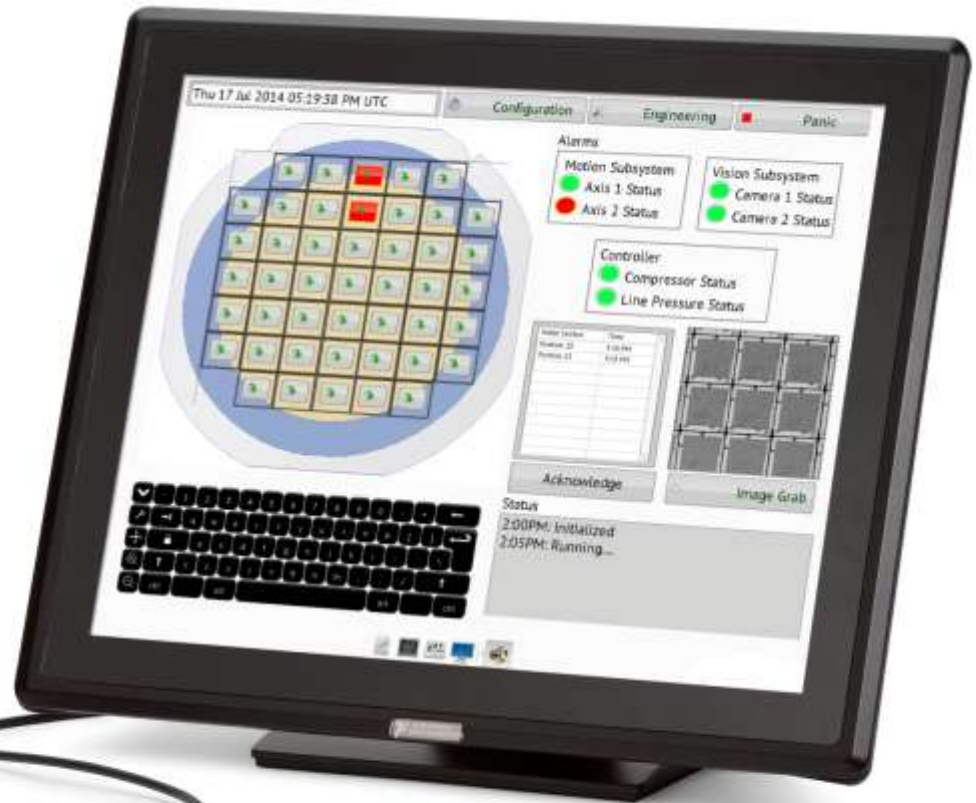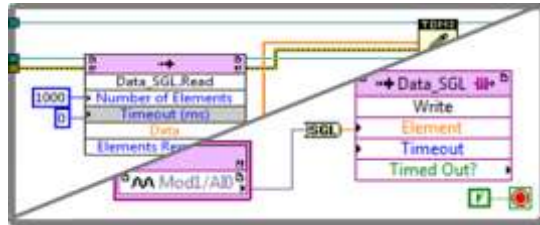


Growing LabVIEW Ecosystem

LabVIEWTools Network

3 million add-ons downloaded
Over 250 add-ons

User Community

Over 10,000 certified users
Over 800 Alliance Partners
Over 80 registered user groups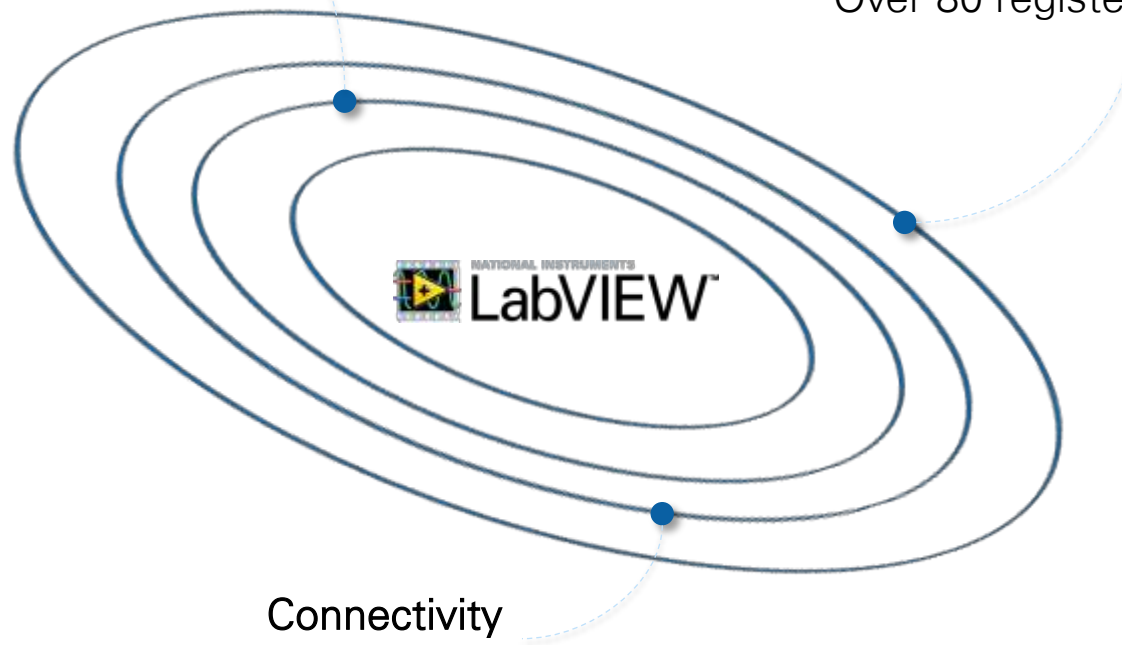

Connectivity

Over 10,000 instrument drivers
Any bus, any protocol, any platform
Integration with third-party languages


LabSocket by Bergmans Mechatronics


New on the LabVIEW Tools Network

- Automatically creates interactive HTML5 interfaces from LabVIEW front panels
- Access on any modern mobile or desktop browser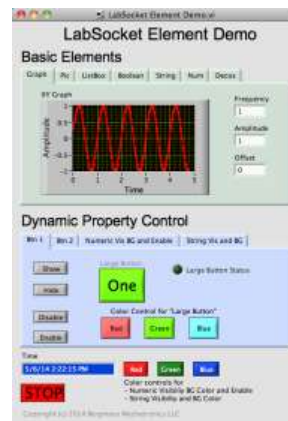

LabVIEW

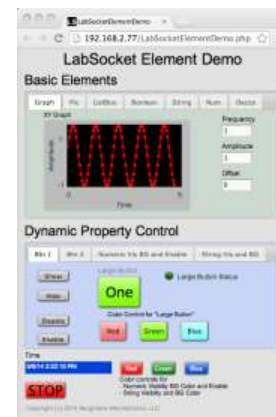Web Browser

Download evaluation at ni.com/labviewtools/labsocket

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